

Bee Stinger

*A Gastight Syringe for the
BASi Bee Syringe Pump*

MD-00001

Laboratory syringe for research purposes. Not for human use.

Important! Before using this syringe remove the plunger from the barrel. Dip plunger into deionized or distilled water and re-insert into glass barrel. Move the plunger up and down five times. Movement of the syringe plunger in a dry barrel may cause excessive wear on the teflon seal, reducing the lifetime of the seal and causing leaks.

Cleaning: Deionized or distilled water, acetone or methyl alcohol are all recommended cleaning solvents for this syringe. Avoid detergents, phosphates and alkalines whenever possible. Syringes may be soaked overnight in a mild solvent or immersed in an ultrasonic cleaner for a few hours if necessary.

765-463-4527
fax 765-497-1102
bas@bioanalytical.com
www.bioanalytical.com



2701 Kent Avenue
West Lafayette, IN
47906 USA